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Types of BSC

Class	Min Inflow Velocity (fpm)	Recirc. Air	Exhaust Air	Contaminated Plenum Surrounded by	Exhaust Alternatives	Biosafety Level
I	75	0%	100%	Outside air (Lab room)	Inside room / Hard Duct	1,2,3
II A1	75	70%	30%	Outside air (Lab room)	Inside room / Thimble Duct	1,2,3
II A2	100	70%	30%	Negative pressure	Inside room / Thimble Duct	1,2,3
II B1	100	30%	70%	Negative pressure	Hard duct only	1,2,3
II B2	100	0%	100%	Negative pressure	Hard duct only	1,2,3
III	Closed P>0.5"WG	0%	100%	Negative pressure	Inside room / Hard Duct	1,2,3,4

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Bio Safety Level (BSL)

	Lethality	Medium	Cure	Example
1	Safe	Liquid	Yes	B.Subtilis
2	Some	Liquid	Some	HIV
3	Serious	Airborne	Some	TBC
4	Extreme	Airborne	None	Ebola

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Table 1. Summary of Biosafety Levels and Infections Agents

Biosafety Level	Infectious Agents	Examples	Practices and Techniques	Safety Equipment (Primary Barriers)	Facilities (Secondary Barriers)
1	<i>Not known to consistently cause disease in healthy adults</i>	Bacillus subtilis, Naegleria gruberi, Infectious canine, hepatitis virus, E.coli	<i>Standard Microbiological practices</i>	None Required	Open benchtop sink required.
2	Associated with human disease. Primary hazards are percutaneous injury, ingestion, mucous membrane exposure	Measles virus, salmonellae, Toxoplasma spp, Hepatitis B Virus.	<i>Level 1 plus:</i> <ul style="list-style-type: none"> Limited access Biohazard warning signs "Sharps" precautions Biosafety manual defining any needed waste decontamination or medical surveillance policies Respiratory protection as required 	<i>Primary barriers:</i> Class I or II BSCs or other physical containment devices used for all manipulations of agents that cause splashes or aerosols of infectious materials. <i>Personal protective equipment:</i> lab coats, gloves, face protection as needed	<i>Level 1 plus:</i> <ul style="list-style-type: none"> Autoclave available
3	Indigenous or exotic agents with potential for aerosol transmission, disease may have serious or lethal consequences	M. Tuberculosis, St. louis encephalitis virus, Coxiella Burnetii, Bacillus anthracis (production level)	<i>Level 2 plus:</i> <ul style="list-style-type: none"> Controlled access Decontamination of waste Decontamination of lab clothing Baseline serum samples of lab personnel 	<i>Primary barriers:</i> Class I or II BSCs or other physical containment devices used for all open manipulations of agents. <i>Personal protective equipment:</i> lab clothing, gloves, respiratory protection as needed	<i>Level 2 plus:</i> <ul style="list-style-type: none"> Physical separation from access corridors Self-closing, double door access Exhaust air not recirculated Negative airflow into laboratory
4	Dangerous / exotic agents that pose a high risk of life threatening disease. Aerosol-transmitted lab infections or related agents with unknown risk of transmission	Ebola zaire, Sin Nombre Virus, Rift Valley Fever	<i>Level 3 plus:</i> <ul style="list-style-type: none"> Clothing change before entry into lab Shower on exit All material decontaminated on exit from lab 	<i>Primary barriers:</i> All procedures conducted in Class III or I or II BSCs in combination with full-body, air-supplied, positive-pressure personnel suit	<i>Level 3 plus:</i> <ul style="list-style-type: none"> Separate building or isolated zone Dedicated supply/exhaust vacuum, and decontamination systems Additional requirements as outlined in Biosafety in Microbiological and Biomedical Laboratories

Source: **Biosafety in Microbiology and Biomedical Laboratories, U.S.** Department of Health and Human Services, HHS publication (CDC) 99-8395. 4th Edition. April 1999. p 52-53.

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Table 2. Biological Safety Cabinet Classes

Source: **Laboratory Biosafety Manual**. Second Edition (revised). Interim guidelines. World Health Organization. p32.

Class	Inflow Velocity (m/s)	Recycle Air (%)	Exhaust Air (%)	Control Plenum Surrounded by	Exhaust Alternatives	Biosafety Level
I	US:0.38 EN:0.70	0	100	Outside Air	Inside room / Hard duct	1, 2 & 3
II Type A1	US:0.38 EN:0.40	70	30	Outside Air	Inside room / Thimble duct	1, 2 & 3
II Type A2	US:0.50 EN:0.40	70	30	Negative Plenum	Inside room / Thimble duct	1, 2 & 3
II Type B1	US:0.50 EN:0.40	30	70	Negative Plenum	Hard duct only	1, 2 & 3
II Type B2	US:0.50 EN:NA	0	100	Negative Plenum	Hard duct only	1, 2 & 3
III	Closed: * >0.5"WC	0	100	Negative Plenum	Inside room / Hard duct	1, 2, 3 & 4

* Pressure differential between chamber and environment.

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
Table 3. *Biological Safety Cabinet Selection*

Type of protection	BSC Selection
Personnel Protection, microorganisms in Risk Groups 1-3	Class I, Class II, Class III
Personnel Protection, microorganisms in Risk Groups 4, glove box laboratory	Class III
Personnel Protection, microorganisms in Risk Groups 4, suit laboratory	Class I, Class II
Product Protection	Class II, Class III only if laminar flow included
Volatile radionuclide / chemical protection, minute amounts	Class II Type B1, Class II Type A2 vented to the outside
Volatile radionuclide / chemical protection	Class I, Class II Type B2, Class III

Source: **Laboratory Biosafety Manual**. Second Edition (revised). Interim guidelines. World Health Organization. p29.

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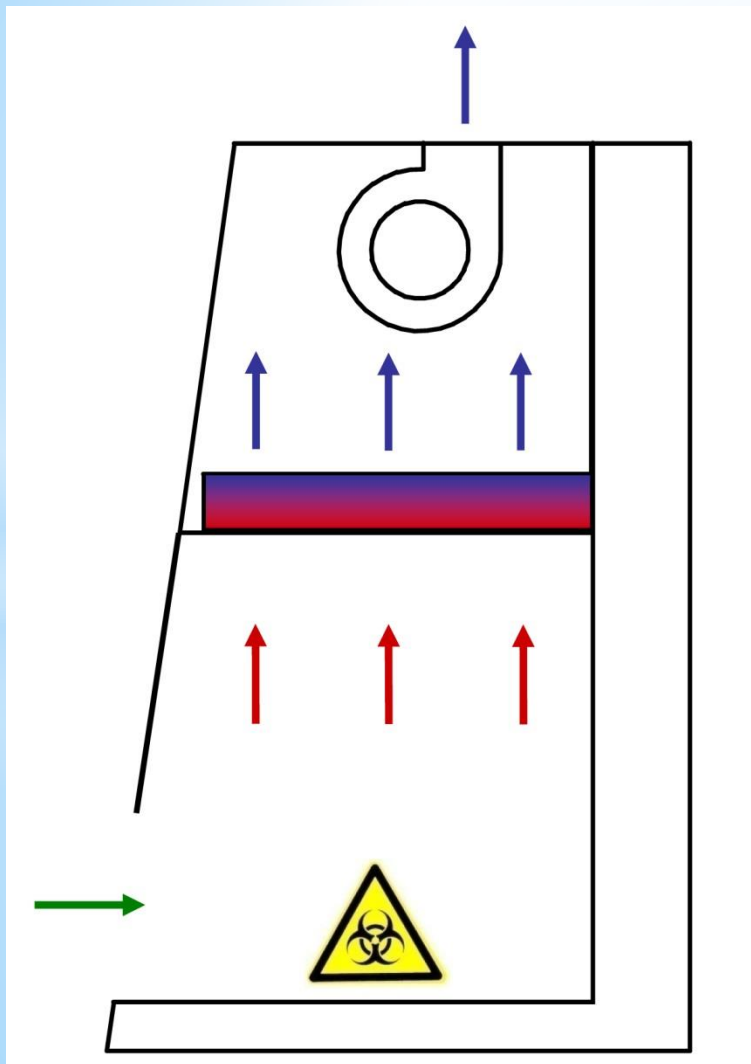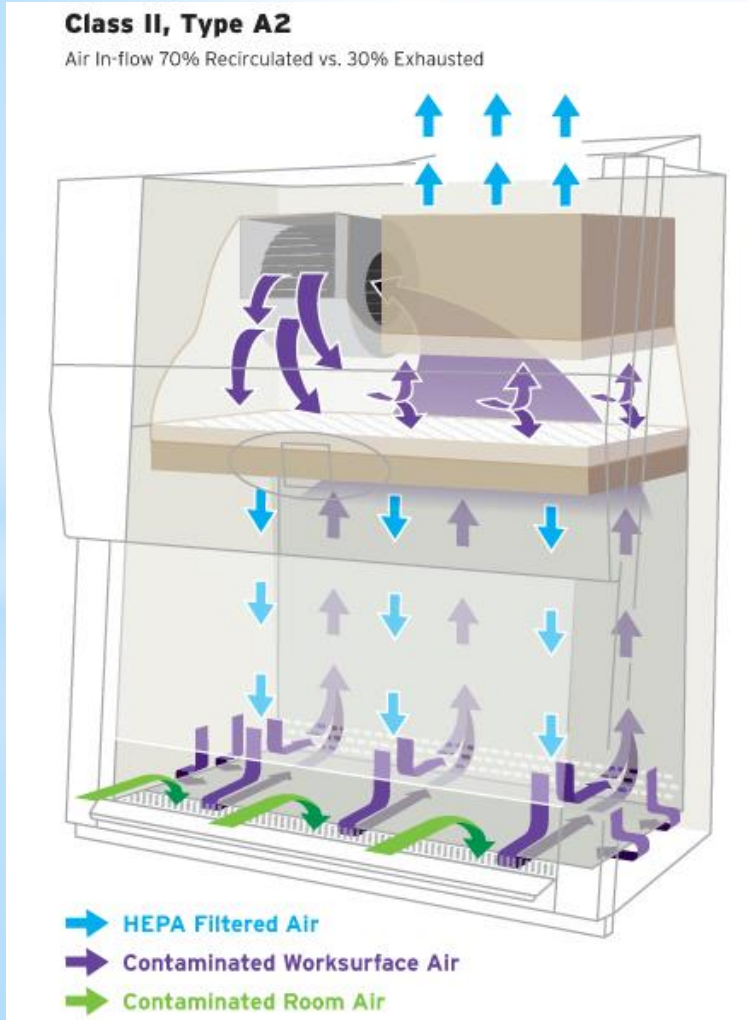
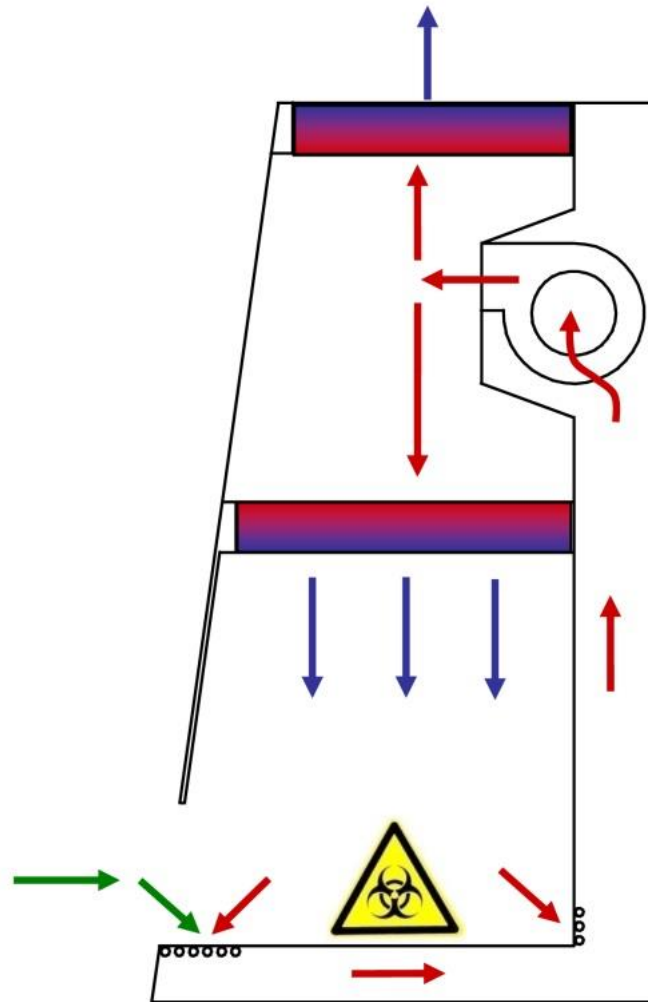
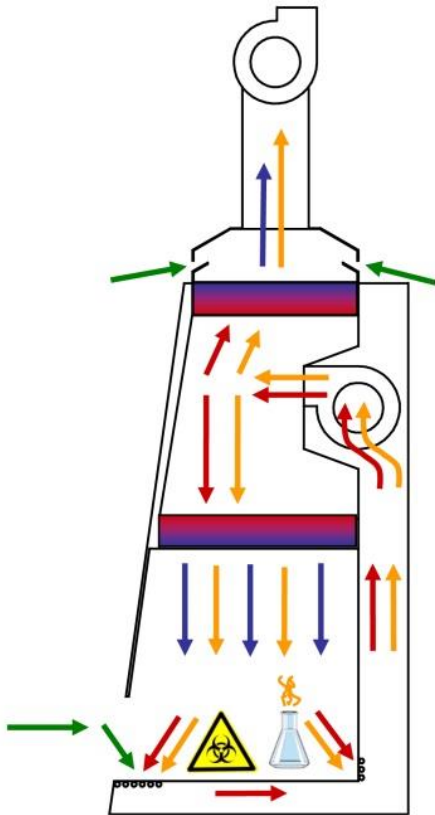
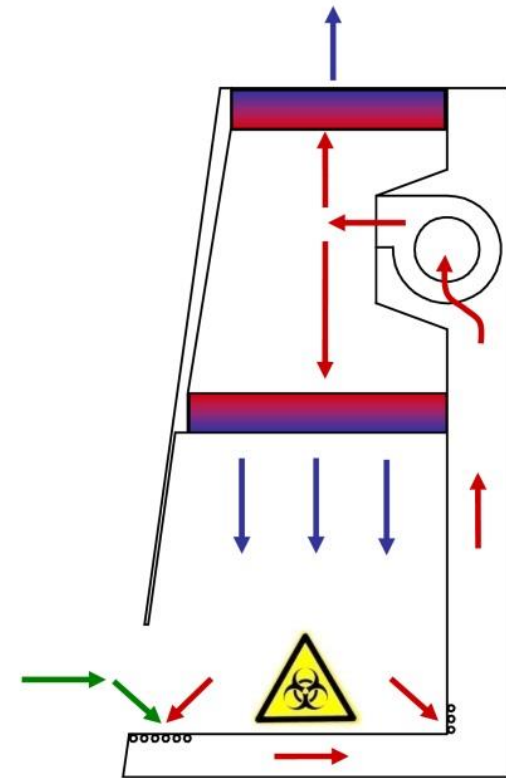
Class II Type A2 with Thimble Ducting

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